

When not in use, keep the accessory in an area (e.g., cabinet or storage case) that is free of excessive dust. If the mirrors do become dusty, use a jet of clean air or inert gas to remove the dust. Do not become concerned with a small amount of dust on any mirror. Such a condition will, generally, have little or no effect on the performance of your attachment.

Avoid touching the mirror surfaces or spilling chemicals on them. Fingerprints and residual chemicals can, in many cases, adversely affect the performance of your attachment.

Avoid contacting the mirror surfaces with metal objects (e.g., wrenches, screwdrivers, spatulas). These may cause scratches or removal of material that could, again, affect the performance.

Should any condition develop with your Harrick accessory that you feel requires contact cleaning and/or refinishing of a mirror, please call our Customer Service Department prior to taking any action. We will be glad to help you.